Finnpipette[®]

Flaschenaufsatz-Dispenser
Bottle-top Dispenser

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Instruction Manual

Please read before use!

Thermo Fisher Scientific Oy
Ratastie 2, P.O.Box 100
FI-01621 Vantaa
Finland

info.pipettes@thermofisher.com
www.thermo.com/finnpipette

1508040

Inhalt

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Contents

	Page
1. Warning and safety Instruction	4
2. Individual components	8
3. Functional Principle	9
4. Prohibited use	10
5. Restrictions on use	11
6. Preparing the unit	12
6.1 Securing the suction tube	12
6.2 Securing the ejection unit	14
6.3 Screwing the unit onto the bottle	15
7. Dispensing	16
7.1 Closure cap removal	17
7.2 Venting the unit	18
7.3 Volume configuration	19
7.4 Dispensing	20
7.5 After use, locking of dispensing mechanism	21
8. Cleaning	22
8.1 Service/ Cleaning	23
8.2 Intensive cleaning	28

9.	<i>Sterilisation</i>	30	9.	<i>Sterilization</i>	30
9.1	<i>Vorbereitung</i>	30	9.1	<i>Preparation</i>	30
9.2	<i>Dosierkolben ausbauen</i>	30	9.2	<i>Removal the piston</i>	30
9.3	<i>Ansaug-/Ausstoßventil lockern</i>	32	9.3	<i>Loosening the suction/ejection valve</i>	32
9.4	<i>Sterilisation durchführen</i>	33	9.4	<i>Sterilization</i>	33
9.5	<i>Dosierkolben wieder einbauen</i>	34	9.5	<i>Reassembly of piston</i>	34
9.6	<i>Ansaug-/Ausstoßventil festdrehen</i>	35	9.6	<i>Tightening the suction/ejection valve</i>	35
9.7	<i>Gerät zusammenbauen</i>	35	9.7	<i>Reassembly of the dispenser</i>	35
10.	<i>Ventile auswechseln</i>	36	10.	<i>Valve replacement</i>	36
10.1	<i>Ansaugventil auswechseln</i>	36	10.1	<i>Suction valve replacement</i>	36
10.2	<i>Ausstoßventil auswechseln</i>	38	10.2	<i>Ejection valve replacement</i>	38
11.	<i>Reparatur beim Hersteller</i>	39	11.	<i>Repair by manufacturer</i>	39
12.	<i>Volumenprüfung</i>	40	12.	<i>Volumetric inspection</i>	40
13.	<i>Problem - Lösungen</i>	42	13.	<i>Problem solutions</i>	44
14.	<i>Technische Daten/ Materialien</i>	46	14	<i>Technical data/ Materials</i>	46
15.	<i>Zubehör und Ersatzteile</i>	49	15.	<i>Accessories and spare parts</i>	49

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1. Warning and safety instructions

Attention!

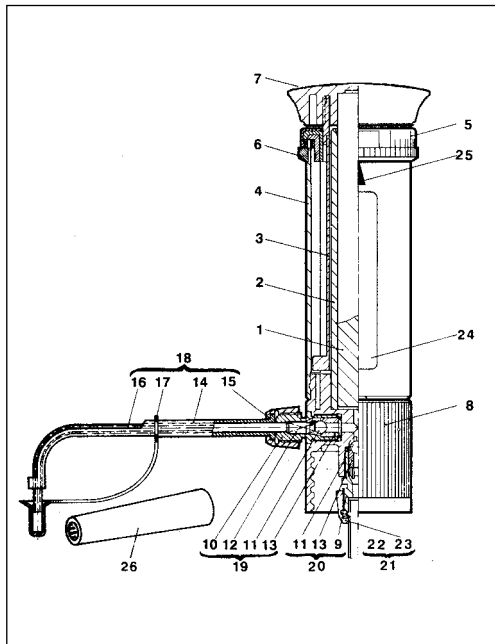
Before use please read the safety instructions and observe them carefully. Before use every member of staff must be familiar with these safety guidelines and they must be kept available at all times. This instruction manual cannot describe all possible safety hazards. It is every user's responsibility to observe all health and safety precautions and to determine and observe restrictions before use of the unit.

- Observe all general safety practices and guidelines for laboratories, such as the use of protective clothing, goggles and gloves when handling hazardous chemicals.

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3. Functional principle

- This instrument is a bottle top dispenser with
digital volume selection for the dispensing of
liquids. The basic model plus the various adap-
ters (see accessories) enable its use with com-
monly used reagent bottles.
- When the piston is raised the ball of the suction
valve is lifted. Liquid is aspirated through the suc-
tion tube from the reservoir bottle into the cylinder.
After finishing the suction phase the suction valve
closes.
- Pushing down the piston opens the ejection valve
and dispenses the volume which was selected on
the volume scale.

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To select the required volume the piston must be in the down position. Turn the adjusting ring in either direction, so that the required volume mark is set above the adjustment arrow.

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4. Prohibited use

The unit should not be used with the following:

- hydrofluoric acid (HF), it attacks the glass.
- liquids that attack Halar (ECTFE), FEP, PTFE, borosilicate glass 3.3, Al_2O_3 or Hastelloy.
- suspensions (e.g. of activated carbon solvent), because solid particles might block the valves.
- easily-combustible media (note the flashpoint).
- use in potentially-explosive environments.

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5. Restrictions on use

- concentrated saline solutions, highly concentrated alkaline solutions may cause the piston to stick or move stiffly.
- the dispenser is suitable for operating temperatures between 4°C and 50°C
- The unit should only be used in areas where it can be assured that no hazardous potentially-explosive atmospheres occur
- Do not deposit in potentially-explosive areas.

Attention - important hint!

Never use force if the piston sticks or becomes stiff. Stop use immediately and carry out a thorough cleaning (see point 8.2).

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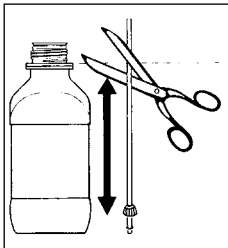
6. Preparing the unit

Attention, please note before assembly!

- For safe and trouble free operation only use the genuine manufacturer's accessories and parts.
- Never use damaged or other than genuine manufacturer's parts.

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
6.1 Securing the suction tube

- Hold the suction tube beside the storage bottle. The tube end with the union nut should face towards the base of the bottle.
- Cut the tube at a slant level to the lower edge of the locking ring on the storage bottle.

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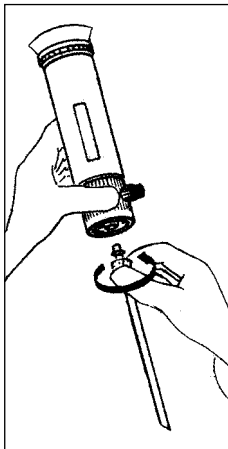
- Push the suction tube on up to the stop.
- Then tighten the union nut onto the valve mantle.

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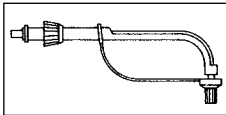
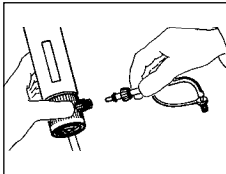
- Put the ejector unit on up to the stop.

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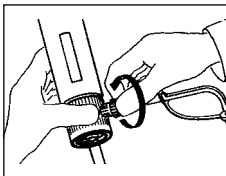- Tighten the union nut firmly and ensure that it is securely fitted.

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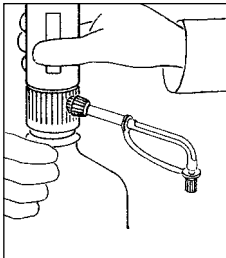
Attention!

- Ensure that the ejection unit is firmly attached
- Only use genuine manufacturer's ejection unit (see accessories and spare parts list).
- Never use deformed or damaged ejection units.

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6.3 Screwing the unit onto the bottle

- Screw the dispenser onto the reservoir bottle, either directly or with one of the enclosed adapters, depending on the bottle's thread size.

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Warning!

- Observe all safety instructions, usage prohibitions and restrictions on use.
- Wear protective clothing, safety glasses and protective gloves.

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7. Dispensing

Attention!

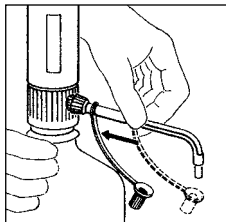
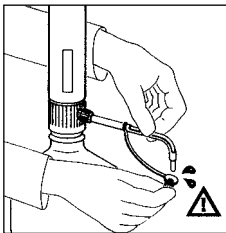
- Follow safety instructions (see point 1).
- Usage prohibitions and restrictions of use must be observed (see point 4 and 5).
- Dispense only when the ejection unit is turned away from the user and any other person.
- Never press down the piston when the closure cap is in fitted to the tube.

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7.1 Closure cap removal

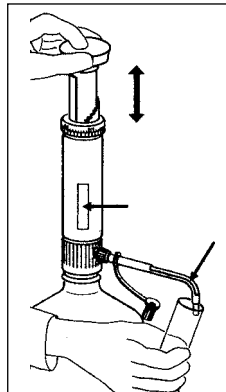
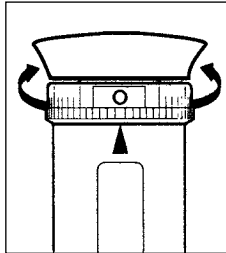
- Position a collecting vessel under the ejection unit.
- Remove closure cap as shown in the drawing.

Caution: liquid might drip off!

- Slide the closure cap's ring to the direction of the device's body as shown in the drawing.

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7.2 Venting the unit

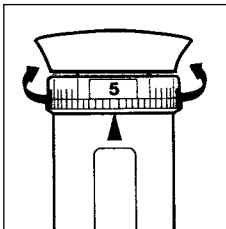
- Turning the scale ring away from the zero position will set the unit for operation.
- Position a collecting vessel under the ejection unit.
- Check whether closure cap is opened. (precautionary)
- Eject air bubbles by pulling the piston a short distance and firmly pressing it down.
- Repeat this operation a few times until no more air bubbles are visible in the observation window (see arrows).
- Start operation (small volume) until no more air bubbles are visible in the ejection tube unit (see arrows).

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- If larger air bubbles can not be ejected, a valve or the suction tube is not fitting tight (see problem solutions point 13).

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7.3 Volume configuration

- By turning the adjusting ring (in either direction) set the desired volume mark directly above the arrow.

Achtung!

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Attention!

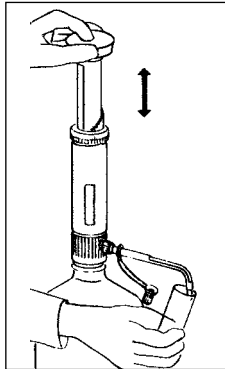
Do this only when the piston is in down position

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7.4 Dispensing

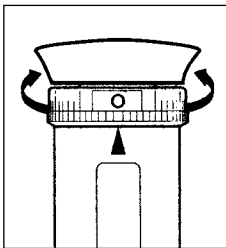
Attention!

Never use force when dispensing. If the piston becomes stiff the unit must be cleaned immediately (see point 8).

- Hold the collecting vessel under the ejection unit.
- Ensure that closure cap is removed.
- Gently and steadily pull up the piston to the top of its path.
- Push the piston down gently and steadily.

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7.5 After use, locking of dispensing mechanism

- After use the piston must always be left in the down position.
- Turning the volume scale to zero position locks the mechanism.

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Attention!

The volume scale can only be turned to the zero position when the piston is in the down position.

- Place the closure cap on the ejection tube's opening.
- After use it is recommended to perform maintenance and cleaning (see point 8.1)

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8. Cleaning

In order to ensure a long life for your dispenser, it should be cleaned regularly.

It is essential to clean it in the following cases:

- piston becoming stiff (see point 8.1)*
- after use with liquids of which the use is restricted (see point 5)*
- before changing reagent*
- before sterilization*
- before replacement of suction or ejection valves*
- before storage*

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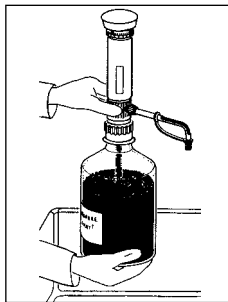
Attention!

- When cleaning always wear protective clothing, goggles and gloves.
- Follow safety instructions (see point 1).
- Suction tube, ejection unit, cylinder and valves contain liquid. Ensure, that they are kept away from your body or other persons.

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
8.1 Service/ Cleaning


- After use set the scale ring to position "0" and close the cap (see point 7.5).
- Place the unit and bottle into a suitable cleaning container

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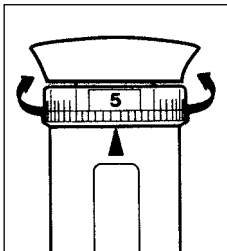- Set the dispenser to maximum volume at the arrow (adjusting mark) through turning the volume scale (left/right).

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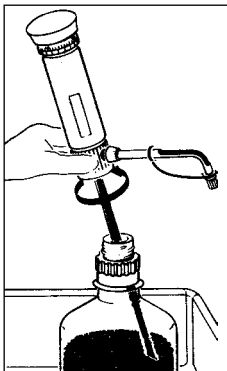- Unscrew the dispenser from the bottle. Ensure that protective gloves are worn.

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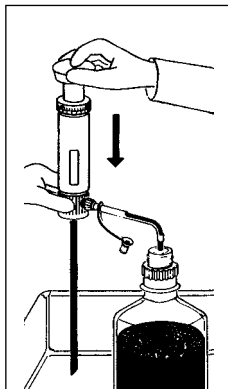
Attention!

Liquid may drip from the suction tube.

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
- Hold the ejection unit above the reservoir bottle.
- Pull off the closure cap and slide it backwards.
- Move the piston up and down to eject the remaining liquid.

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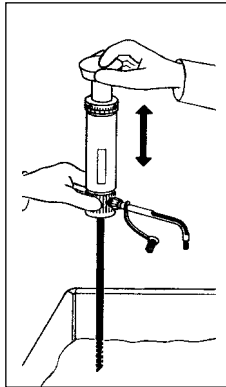
- Hold the suction tube into a suitable cleaning solution and rinse the dispenser completely by repeated dispensing.

Attention: Liquid may splash!
Always dispense away from your body!

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
- Rinse the unit with suitable solvent by repetitive dispensing.
- Remove the suction tube from the solvent and clear residual solvent from the dispenser by repeated dispensing.

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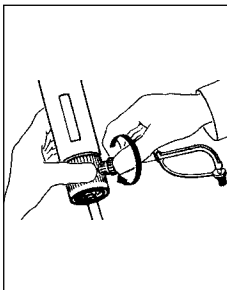- Dismantle the dispenser and clean all following parts separately after dismantling the dispenser as described following:
- Unscrew the ejection unit from the dispenser.

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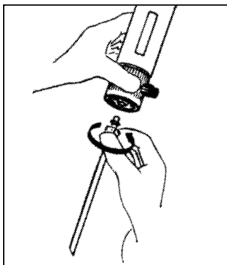
- Remove the union nut on the suction tube.
- Remove the suction tube from the dispenser.

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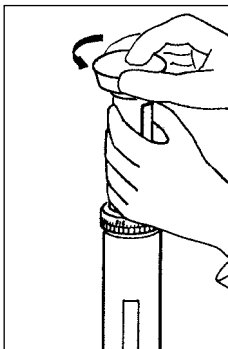
8.2 Intensive cleaning

- Clean the unit as described in point 8.1.
- Remove the piston as follows:
 - Set the maximum volume at the arrow by turning the adjusting ring.
 - Pull the piston right to the stop.
 - Hold the step gauge and unscrew the knob/piston.
 - Pull the piston out of the cylinder.

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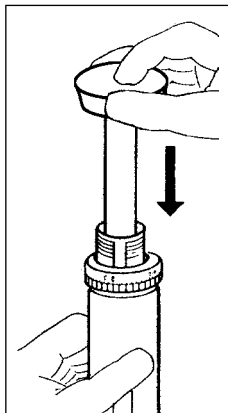- Clean the cylinder and the ceramic piston with a brush and detergent. Afterwards rinse these parts with an appropriate solvent.
- Reassemble the piston: Keep the unit vertical and insert the piston carefully into the cylinder, with a **smooth rotating** movement. Screw the knob (with the piston) onto the step gauge.

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Attention!

Take care that the piston is inserted straight and no force is used.

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9. Sterilization

The dispenser is able to withstand autoclaving up to 121 °C, 2 bar for a period of time acc. to DIN te = 20 min.

9.1 Preparation

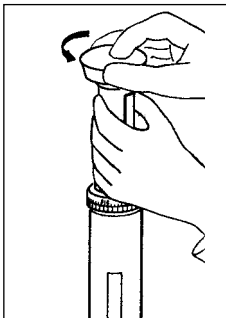
- Before sterilization clean the unit as described in point 8.1.

9.2 Removal of piston

Hint!

In order to achieve a better cleaning effect, remove the piston from the cylinder

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

- Set the maximum volume at the arrow, by turning the adjusting ring.
- Pull the piston right to the stop.
- Hold the step gauge and unscrew the knob/piston.
- Pull the piston out of the cylinder .

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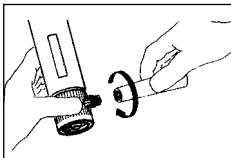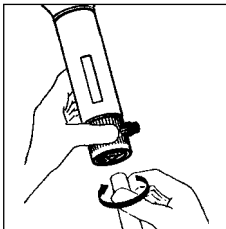
9.3 Loosening the suction/ ejection valve

- Loosen the suction valve with the assembly wrench by turning it counterclockwise a half a revolution.
- Loosen the ejection valve with the assembly wrench by turning it counterclockwise a half a revolution.

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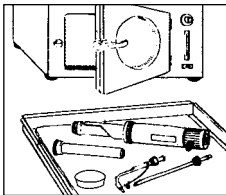9.4 Sterilization

Place all parts in the autoclave (up to 121 °C, 2 bar, for a period of time acc. to DIN te = 20 min.)

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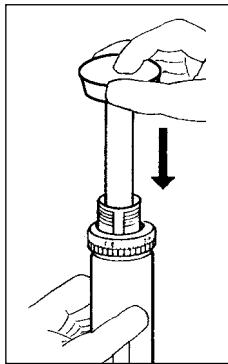
Attention!

- Let the unit cool down to room temperature, before reassembling.
- Check parts for damage and if necessary, exchange them.

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9.5 Reassembly of piston

- Keep the unit vertically and insert the piston **carefully** into the glass cylinder with a **smooth rotating** movement.

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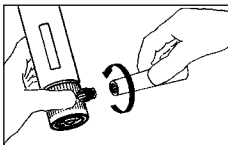
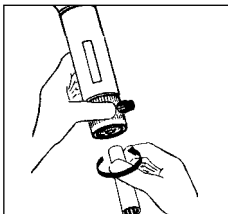
9.6 Tightening the suction/ ejection valve

- Tighten the suction valve firmly with the assembly wrench by turning it in a clockwise direction.
- Tighten the ejection valve firmly with the assembly wrench by turning it in a clockwise direction.

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9.7 Reassembly of the dispenser

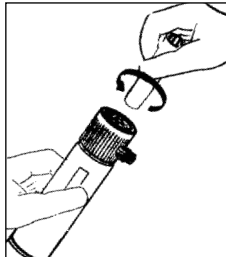
- Reassemble the unit (see point 6) and check correct operation (see point 7). In case of problems see the trouble shooting guide (p. 13).

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
10. Valve replacement

Leaking or damaged suction/ejection valves require immediate exchange (see accessories and spare parts point 15).

10.1 Suction valve replacement

- *Clean the unit (see point 8).*
- *Unscrew the suction tube.*
- *Hold the unit upside down, otherwise the valve ball might fall out, when changing the valve.*
- *Unscrew the suction valve with the assembly wrench and remove the valve ball.*

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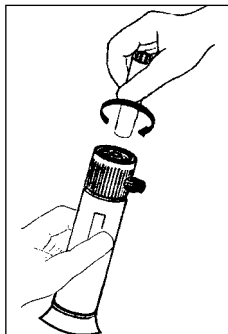
- Insert the ball again and screw on the new suction valve, tightening with the assembly wrench.

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- Put in the suction tube right to the stop and tighten the union nut on the valve (see point 6.1).

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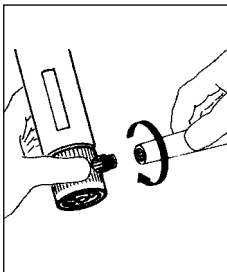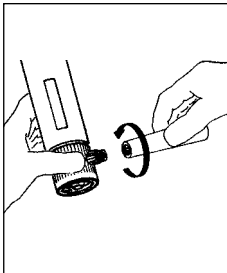
10.2 Ejection valve replacement

- Realise cleaning (see point 8).
- Unscrew the ejection unit (see point 8).
- Screw out the ejection valve with the assembly wrench and remove.

- Screw in the new ejection valve and tighten firmly with the assembly wrench.

- Push on the ejection unit up to the stop and tighten the union nut firmly (see point 6.2). Check to ensure a ensure a secure fitting!

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11. Repair by manufacturer

- A defective unit should be cleaned carefully prior to sending for repair (see chapter 8).
- Units will only be repaired if they are cleaned by the user and are no longer contaminated with any substances which could pose a risk to health.
- The clearance certificate should be filled out prior to sending in the unit (see point 16).
- Not cleaned units will be returned to the sender without having been repaired.
- Describe the malfunction exactly.
- Indicate the dispensed liquid. (Important!)
- The warranty is invalid in case unit has been maintained other than described in this manual (e.g. use of force).

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12. Volumetric inspection

The terms for accuracy (R%) and the coefficient of variation (CV%), indicated in point 14, can be checked as follows:

1. Measuring procedure

(see e.g. DIN EN ISO 8655-6):

- Set the nominal volume
 - Dispense with de-ionized water
 - Determine the dispensed quantity on an analytical balance
 - Convert the measured weight to volume
- Note:** Temperature, air pressure and material need to be considered

2. Repetitive measuring

- In order to obtain reliable measures for accuracy and coefficient of variation, the procedure described under 1. should be repeated at least 5 times.

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3. Calculation of accuracy and coefficient of variation

- Accuracy (R%) and coefficient of variation (CV%) have to be calculated in accordance to the formulas of statistical quality control.

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13. Problem solutions

<i>Symptom</i>	<i>Possible cause</i>	<i>Solution</i>
<i>Piston is stiff or does not move at all</i>	<ul style="list-style-type: none">- Closure cap not open- Crystallisation at the piston or the valves	<ul style="list-style-type: none">- Open the closure cap (see point 7.1) unit- Stop dispensing and clean the thoroughly (see point 8.2). If necessary, soak it for 24 hours in a bath (e.g. with 20 % reasol cleaning solvent).
<i>Suction or dispensing is impossible</i>	<ul style="list-style-type: none">- Volume scale set on "0" position- Suction/ejection valve sticking	<ul style="list-style-type: none">- Turn the adjusting ring to set the required volume (see point 7.2 and 7.3).- Clean the unit (see point 8.1)- Shake the unit or tap the valve block against the palm of the hand.- If necessary, change the valves (see point 10).
<i>Liquid is leaking from the gap between the valve block and the valves.</i>	<ul style="list-style-type: none">- Valves loose or damaged	<ul style="list-style-type: none">- Tighten valves with the assembly wrench Note: After sterilization (see point 9), suction/ejection valves need to be tightened.

<i>Liquid is leaking between the ejection unit and the valve</i>	<ul style="list-style-type: none"> - Ejection unit is loose or damaged - Wrong ejection unit is used 	<ul style="list-style-type: none"> - Mount the original ejection unit correctly (see point 6.2). - Use only the original ejection unit.
<i>Unit is aspirates in air</i>	<ul style="list-style-type: none"> - Unit is not primed - Suction tube is loose - Suction tube or union nut damaged 	<ul style="list-style-type: none"> - Prime the unit (see point 7.2). - Tighten the suction tube (see point 6.1). - Replace original suction tube (see accessories spare parts point 15).
<i>Delivered volume is incorrect</i>	<ul style="list-style-type: none"> - Ejection unit loose or damaged - Suction tube loose or damaged - Suction/ejection valves loose or damaged 	<ul style="list-style-type: none"> - Install a geneime ejection unit correctly (see point 6.2). - Tighten the suction tube (see point 6.1). - Tighten the valves with the assembly wrench. - Clean the unit (see point 8.1). - If necessary, change the valves (see point 10).

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Dispenser Volumen ml		Einstell- volumen ml	R* %	CV* %
0,2	- 1	0,05	* 0,5	* 0,2
0,4	- 2	0,05	* 0,5	* 0,1
1	- 5	0,10	* 0,5	* 0,1
2	- 10	0,25	* 0,5	* 0,1
5	- 30	0,50	* 0,5	* 0,1
10	- 60	1,00	* 0,5	* 0,1

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* The given values for accuracy and coefficient of variation refer to the maximum volume at
 - equal temperature (20 °C) of the unit, surrounding and liquid (deionized water)
 - regular, non-jerky operation

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The dispenser is conformity certified acc. to the prescription of the German "Eichordnung" (Enclosure 12).

Area of temperatures for use of the unit and the liquid between 4 °C and 50 °C.

Material

Material

<i>Kolben</i>	<i>Zylinder</i>	<i>Ventil-</i>	<i>Ventil-</i>	<i>Ventil-</i>	<i>Ventil-</i>
<i>piston</i>	<i>cylinder</i>	<i>block</i>	<i>kugel</i>	<i>sitz</i>	<i>feder</i>
		<i>valve</i>	<i>valve</i>	<i>valve</i>	<i>valve</i>
		<i>block</i>	<i>ball</i>	<i>seat</i>	<i>spring</i>

<i>höchstreine</i>	<i>Borosilikat-</i>	<i>Halar</i>	<i>Borosilikat-</i>	<i>Borosilikat-</i>	<i>Hastelloy</i>
<i>Al₂O₃-Keramik</i>	<i>Glas 3.3</i>	<i>(ECTFE)</i>	<i>Glas 3.3</i>	<i>Glas 3.3</i>	
<i>(99,7%)</i>					

<i>highly pure</i>	<i>Borosilicate</i>	<i>Halar</i>	<i>Borosilicate</i>	<i>Borosilicate</i>	<i>Hastelloy</i>
<i>Al₂O₃-ceramic</i>	<i>glass 3.3</i>	<i>(ECTFE)</i>	<i>glass 3.3</i>	<i>glass 3.3</i>	
<i>(99,7%)</i>					

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CAUTION!

The Finnpipette is designed to allow easy in-lab service. If you would prefer to have us or your local representative service your pipette, please make sure that the pipette has been decontaminated before you send it to us.

Please note that the postal authorities in your country may prohibit or restrict the shipment of contaminated material by mail.