Thermo Scientific S1 Pipet Filler

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Disposal of the Pipette

Follow the guidelines below for the disposal of the S1 Pipet Filler.

Warning Decontaminate the pipette before disposal.

Follow laboratory and country-specific procedures for biohazardous or radioactive waste disposal.

Warning The used lithium (Li-ion) battery is regulated waste and must be disposed of according to local regulations.

Dispose of the instrument according to the legislation stipulated by the local authorities concerning take-back of electronic equipment and waste.

The procedures vary by country.

Pollution degree: 2 according to IEC 60664-1 *

Method of disposal:

- Electronic waste
- Contaminated waste
- (Infectious waste)

Regarding the original packaging and packing materials, use the recycling operators known to you. For more information, contact your local Thermo Fisher Scientific representative.

*) The pollution degree describes the amount of conductive pollution present in the operating environment. Pollution degree 2 assumes that normally only nonconductive pollution, such as dust, occurs with the exception of occasional conductivity caused by condensation.

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Product specifications are subject to change without prior notice. All trademarks are the property of Thermo Fisher Scientific Inc. and its subsidiaries.

See the latest version at **www.thermofisher.com**

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
Thermo Scientific S1 Pipet Filler

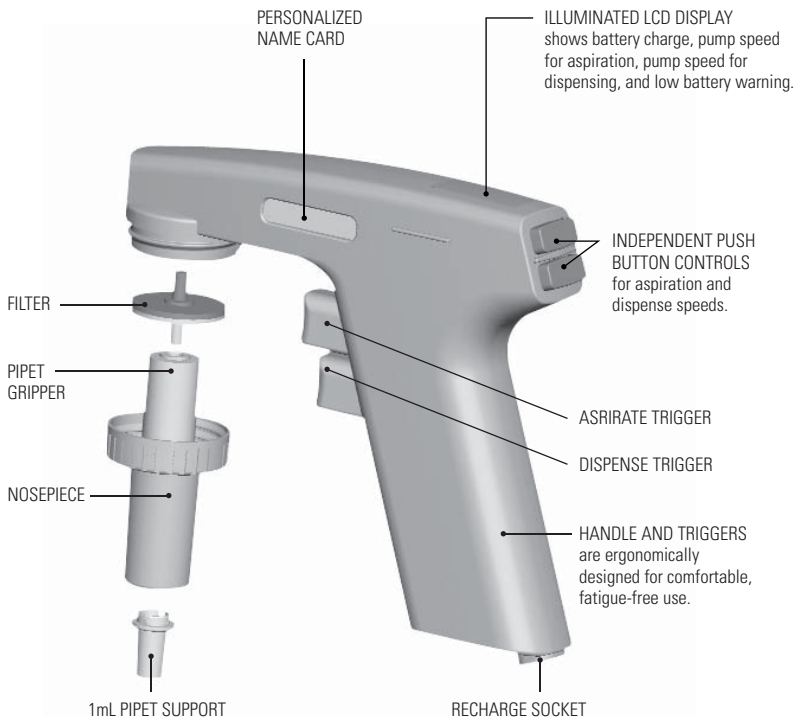Pipettor Overview

MANY USEFUL ACCESSORIES are included to provide for your specific applications. See inside for details.

Thermo Scientific S1 Pipet Filler: Cordless Pipettor With Advanced Solid State Technology

Congratulations!

You have purchased a lightweight, compact, cordless Thermo Scientific S1 Pipet Filler. Designed for use with graduated and volumetric glass and plastic pipets, S1 Pipet Filler offers simple, efficient pipetting performance and maximum pipetting comfort.

Operating on a heavy-duty, rechargeable lithium-ion battery, this pipettor can operate approximately 3 to 4 times longer between charges than typical units using nickel cadmium batteries.

Before You Begin

Included in this S1 Pipet Filler package, you should find the following:

- Thermo Scientific S1 Pipet Filler
- Power Supply (for charging battery or for use with AC current)
- Wall-Mounted Holder
- Table-Mounted Stand
- Pipet Support to Eliminate Wobble with 1mL Pipets
- Name Cards and Clear Covers for Personalizing Your S1 Pipet Filler
- Certificate of Warranty and Warranty Registration Card
- Power Adapters for International Use

If you are missing any of the above items, please contact your Thermo Scientific representative.

Thermo Scientific S1 Pipet Filler

Charging/Recharging

Warning: Use only the original Thermo Scientific S1 Pipet Filler charger and battery pack.

The pipet filler battery may be discharged when delivered and must be charged before initial use. Insert the country specific power plug adapter into the power supply unit. Plug the power supply into the recharge socket at the base of the pipettor's handle and the other end to an AC wall socket. Allow the unit to charge anywhere from 15 minutes to 2 hours. **To extend the battery life cycle, it's recommended to charge the pipet filler every two months even if it is not used on a daily basis. A typical S1 Pipet Filler battery lasts for approximately 400 charging cycles. Please note that a depleted battery can cause the pipette to malfunction. We recommend that at a minimum you should charge the battery every 3 years.** For added ease and convenience, S1 Pipet Filler features:

- A battery icon on the LCD display that indicates the level of battery charge -- the more bars, the greater the charge.
- Special circuitry that makes it impossible to overcharge the battery and that allows the pipettor to be charged or partially charged at any convenient time.

Using S1 Pipet Filler

To use your S1 Pipet Filler, insert a pipet securely in the holder and follow these simple directions:

To Aspirate

Insert the pipet into your sample and depress the upper trigger.

To Dispense

Insert the pipet into your chosen vessel and depress the lower trigger. (Partially depressing the trigger allows for controlled gravity dispensing.)

To Set The Pump Speed

Buttons on the rear of your S1 Pipet Filler unit allow you to set your aspiration (upper button) and dispense (lower button) speeds. Eight speeds are available for each function – simply push “+” for faster and “-” for slower. Your selection will be displayed on the lower portion of the pipettor's LCD. A zero speed selection on dispense enables gravity dispensing.

Disinfecting And Disassembling

The entire S1 Pipet Filler can be cleaned with most common laboratory disinfectants but must not be submerged in liquid. To prepare the unit for cleaning:

- Remove the nosepiece by turning the housing in a counterclockwise direction.
- Remove the silicone rubber pipet gripper from the nosecone. Pipet gripper and nosecone may be autoclaved.

Air Filter

Aspiration rate slowing may be a sign that the air filter needs replacing. To do so, remove the nosepiece and pipet gripper, pull the filter firmly and wiggle to remove from body.

Wall Attachment

The enclosed S1 Pipet Filler wall mount may be attached to a smooth, clean surface using the adhesive backing (also included). For best results, wipe the wall surface with alcohol, let dry, press wall mount firmly in place and allow to set overnight before using. For rough surfaces, mount with screws.

1mL Pipet Support

To enhance pipetting performance when using 1mL pipets, S1 Pipet Filler features a unique pipet support. Simply snap the accessory into the nosecone housing, insert your pipet, and perform your chosen application with a stable, wobble-free pipet.

Power Requirements

Power supply input	100V AC or 240V AC, 50-60 Hz, 0.2 - 0.6 A
Power supply output	6.0V DC, max 1.2 A Constant Current

EXPLANATION OF SYMBOLS ON POWER SUPPLY

---	=	DC Symbol
	=	For Indoor Use Only
	=	Class 2 Power Supply

Environmental Operating Conditions

For optimum performance, your S1 Pipet Filler should be used under the following range of conditions:

Operating altitude	0-1,524m
Operating temperature	10-35°C (For Indoor Use Only)
Storage temperature	-10 to 40 ° C
Relative humidity	10-95% non-condensing
Mains supply voltage fluctuations (for external power supply)	90V AC or 264V AC
Transient overvoltage (for the power supply)	Installation Category II
Applicable rated pollution degree	Pollution Degree 2 (normally only non-conductive pollution occurs)

Thermo Scientific S1 Pipet Filler

Hazards And Precautions

If the equipment is used in a manner not specified by the manufacturer, the protection provided by the equipment may be impaired.

Warning Ensure that the power transformer is connected to a power receptacle that provides the voltage and current specified in this user manual.

Use of an incompatible power receptacle can cause shock and fire hazard.

Warning Incorrect handling of the battery or pipette may cause personal injury. Do not drop the battery. Do not expose the battery to any kind of mechanical stress or extreme temperatures: above 60°C (45°C when charging) or under 0°C.

Warning The S1 Pipet Filler contains a Li-ion battery and a potential malfunction of the battery (short circuit, mechanical damage, overheating, etc.) may cause explosion. Do not touch a leaking battery directly. Do not use the battery in areas with abundant static electricity, otherwise the safety devices may be damaged, creating the potential for hazardous situations.

Caution The charger's plug must be inserted in to a socket that is easy to reach and from which it can be easily unplugged in case of emergency.

Warning Use of incorrect power supply units may result in fatal injuries and damage to the device.

Technical Information

For technical assistance or to reorder, contact your nearest Thermo Scientific representative. See back cover for details.


S1 Pipet Filler models and Accessories

Item No.	Description	Unit
95*	S1 Pipet Filler	1
9066	Single Pipet Filler Wall Mount Stand	1
9067	Single Pipet Filler Table Mount Stand	1
9069	Single Pipet Filler Wing Stand	1
9065	Silicone Pipet Gripper	4/case
9070	1 ml Pipet Filler Support	1
9068	Power Supply (Universal Input)	1
9057	Hydrophobic Filters, 0.45 µm, Sterile	25/bag
4580560	Hydrophobic filters, 0.45 µm, Non-Sterile	5/bag
9064	Nosepiece (holder, silicone gripper, filter)	1

*Available Colors: 01 = White, 11 = Clear, 21 = Blue, 31 = Red, 41 = Green

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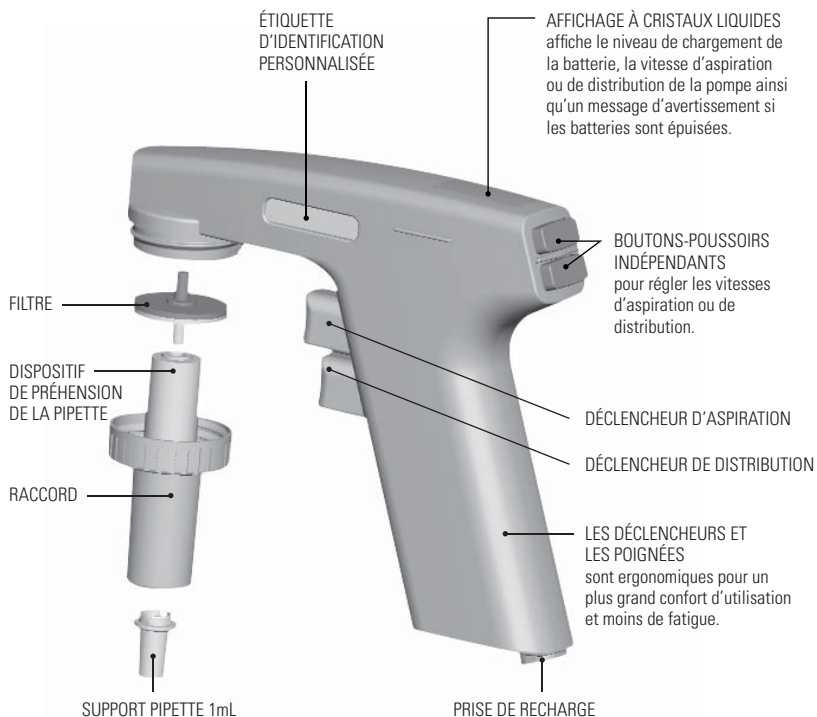Technical Support:

Europe and Asia:
info.pipettes@thermofisher.com

North America:
ushudtechnicalservices@thermofisher.com

Visit www.thermofisher.com for additional product information

Visit www.thermoscientific.com/glp for application notes, videos and to learn good laboratory pipetting practices

thermofisher.com

info.pipettes@thermofisher.com

© 2016 Thermo Fisher Scientific Inc. All rights reserved. All trademarks are the property of Thermo Fisher Scientific Inc. and its subsidiaries. Specifications, terms and pricing are subject to change. Not all products are available in all countries. Please consult your local sales representative for details.

Thermo Fisher Scientific

Ratastie 2, P.O.Box 100
FI-01621 Vantaa
Finland

1508880-03

Thermo
SCIENTIFIC

Part of Thermo Fisher Scientific